



Spike RBD (SARS-CoV-2) : ACE2 Inhibitor Screening Assay Kit

Product Information

Cat.No. Kit-5200

Product Overview

The SARS-CoV-2 Spike:ACE2 Inhibitor Screening Assay Kit is designed for screening and profiling inhibitors of this interaction. This kit comes in a convenient 96-well format, with Hislabeled ACE2, purified SARS-CoV-2 Spike protein, HRP-labeled anti-His antibody, and assay buffer for 100 binding reactions. The key to this kit is the high sensitivity of detection of bound Hislabeled ACE2 by HRP-labeled Anti-His. Only a few simple steps on a microtiter plate are required for the assay. First, SARS-CoV-2 Spike is coated on a 96-well plate. Next, ACE2-His is incubated with SARS-CoV-2 Spike on the plate. Finally, the plate is treated with Anti-His-HRP followed by addition of an HRP substrate to produce chemiluminescence, which then can be measured using a chemiluminescence reader.

Description

The pandemic coronavirus disease 2019 (COVID-19) is caused by Severe Acute Respiratory Syndrome Coronavirus 2 (SARS-CoV-2). As a first step of the viral replication strategy, the virus attaches to the host cell surface before entering the cell. The Spike protein receptor binding domain (RBD) recognizes and attaches to the Angiotensin-Converting Enzyme 2 (ACE2) receptor found on the surface of type I and II pneumocytes, endothelial cells, and ciliated bronchial epithelial cells. Drugs targeting the interaction between the Spike protein of SARSCoV-2 and ACE2 may offer some protection against the viral infection.

Stability

Up to 6 months from date of receipt, when stored as recommended.

Size

96 reactions

Kit Components

SARS-CoV-2 Spike Protein (RBD), mFc Tag, 5 µg, -80°C

ACE2, His-Tag, 5 µg, -80°C

3x Immuno Buffer 1, 50 ml, -20°C



Spike RBD (SARS-CoV-2) : ACE2 Inhibitor Screening Assay Kit

Blocking Buffer 2, 50 ml, +4°C

Anti-His-HRP, 15 µl, -80°C

ELISA ECL substrate A (transparent bottle), 6 ml, Room Temp

ELISA ECL substrate B (brown bottle), 6 ml, Room Temp

96-well white microplate, 1, +4°C

Avoid multiple freeze/thaw cycles!

Materials Required but Not Supplied

PBS (Phosphate buffered saline)

Luminometer or fluorescent microplate reader capable of reading chemiluminescence

Adjustable micropipettor and sterile tips

Assay Protocol

All samples and controls should be tested in duplicate.

Coating the plate with SARS-CoV-2 Spike:

- 1) Thaw SARS-CoV-2 Spike on ice. Upon first thaw, briefly spin tube containing SARS-CoV2 Spike to recover the full contents of the tube. Aliquot into single use aliquots. Immediately store remaining SARS-CoV-2 Spike in aliquots at -80°C. Note: SARS-CoV2 Spike is very sensitive to freeze/thaw cycles. Avoid multiple freeze/thaw cycles.
- 2) Dilute SARS-CoV-2 Spike to 1 µg/ml in PBS.
- 3) Add 50 µl of diluted SARS-CoV-2 Spike solution to each well and incubate overnight at 4°C.
- 4) Dilute 3x Immuno Buffer 1 to 1x Immuno Buffer 1 with water. Dilute only the amount required for the assay; store remaining 3x Immuno Buffer 1 undiluted.
- 5) Decant to remove supernatant. Wash the plate three times with 100 µl 1x Immuno Buffer 1. Tap plate onto clean paper towels to remove liquid.
- 6) Block wells by adding 100 µl of Blocking Buffer 2 to each well. Incubate for 1 hour at room temperature with slow shaking. Remove supernatant as described in step 5.

Step 1:

- 1) Add 20 µl of 1x Immuno Buffer 1 to each well.
- 2) Add 10 µl of inhibitor solution to each well designated "Test Inhibitor". For the "Positive Control"



Spike RBD (SARS-CoV-2) : ACE2 Inhibitor Screening Assay Kit

and “Blank”, add 10 µl of the same solution without inhibitor (inhibitor buffer). Optionally, incubate at room temperature for one hour with slow shaking.

Note: It is recommendable to use PBS to dilute antibodies or other proteins acting as neutralization inhibitors. When using small molecules dissolved in DMSO, final DMSO concentration in the assay should be $\leq 1\%$. Inhibitor buffer should contain the same concentration of DMSO as the test inhibitor.

3) Thaw ACE2-His on ice. Upon first thaw, briefly spin tube containing enzyme to recover full contents of the tube. Aliquot ACE2-His into single use aliquots. Immediately store remaining undiluted enzyme in aliquots at -80°C . Note: ACE2-His is very sensitive to freeze/thaw cycles. Do not re-use thawed aliquots or diluted enzyme.

4) Dilute ACE2-His to 2.5 ng/µl (approximately 30 nM) in 1x Immuno Buffer 1. Keep diluted protein on ice until use. Discard any unused diluted protein after use.

5) Add another 20 µl of 1x Immuno Buffer 1 to the wells designated “Blank”.

Blank Positive Control Test Inhibitor

1x Immuno Buffer 1 40 µl 20 µl 20 µl

Test Inhibitor - - 10 µl

5% DMSO in water (Inhibitor buffer) 10 µl 10 µl -

ACE2-His (2.5 ng/µl) - 20 µl 20 µl

Total 50 µl 50 µl 50 µl

6) Initiate reaction by adding 20 µl of diluted ACE2-His (see Step 1-4) to wells labeled “Positive Control” and “Test Inhibitor”. Incubate at room temperature for one hour with slow shaking.

7) Decant to remove supernatant. Wash the plate 3 times with 100 µl/well 1x Immuno Buffer 1. Tap plate onto clean paper towels to remove liquid.

8) Block wells by adding 100 µl of Blocking Buffer 2 to each well. Incubate for 10 minutes at room temperature. Remove supernatant as in Step 1-7.

Step 2:

1) Dilute Anti-His-HRP 1000-fold with Blocking Buffer 2.

2) Add 100 µl to each well. Incubate for 1 hour at room temperature with slow shaking.

3) Wash plate three times with 1x Immuno Buffer 1. Tap plate onto clean paper towel to remove liquid.



Spike RBD (SARS-CoV-2) : ACE2 Inhibitor Screening Assay Kit

- 4) Block wells by adding 100 µl of Blocking Buffer 2 to each well. Incubate for 10 minutes at room temperature. Decant to remove supernatant. Tap plate onto clean paper towels to remove liquid.
- 5) Just before use, mix 50 µl ELISA ECL Substrate A and 50 µl ELISA ECL Substrate B, then add 100 µl to each well. Discard any unused chemiluminescent reagent after use.
- 6) Immediately read sample in a luminometer or microtiter-plate capable of reading chemiluminescence. "Blank" value is subtracted from all readings.

Reading Chemiluminescence:

Chemiluminescence is the emission of light (luminescence) which results from a chemical reaction. The detection of chemiluminescence requires no wavelength selection because the method used is emission photometry and is not emission spectrophotometry.

To properly read chemiluminescence, make sure the plate reader is set for LUMINESCENCE mode. Typical integration time is 1 second; delay after plate movement is 100 milliseconds. Do not use a filter when measuring light emission. Typical settings for the Synergy 2 BioTek plate reader are: use the "hole" position on the filter wheel; Optics position: Top; Read type: endpoint. Sensitivity may be adjusted based on the luminescence of a control assay without enzyme (typically we set this value as 100).